

BIOBASE®

Portable Refrigerator

LCX-6L/12L/20L

CR-19

User Manual

BIOBASE GROUP

Version 2021.04

Application Method

1. Before using our Portable Refrigerator, carefully check whether the box body is damaged, whether there are holes, whether the buckle is damaged, to ensure that the sealing of the Portable Refrigerator is intact
2. Take out the cold storage agent configured in the box, check whether the cold storage agent is in good condition, and inject water to the marked water injection line;
3. Separate the water-filled cold storage agent, put it into the cold storage or freezer, reserve space appropriately (because the volume of the cold storage agent will expand slightly after freezing), and refrigerate for more than 6-10 hours (the length of time is related to the freezing capacity of the equipment) until the cold storage agent is completely frozen into solid;
4. Put the frozen refrigerant into the Portable Refrigerator;
5. After the temperature in the box reaches the requirement, put the articles in. If there is a partition, please separate the articles from the refrigerant with a partition.
6. Close the lid and check whether the lid is sealed

Attentions

1. After using (for a period of time; when the temperature in the box is higher than 8°C, please replace the completely frozen refrigerant.
 2. After using the Portable Refrigerator, just wipe the box body gently with a dry cloth. Please take good care of the Portable Refrigerator for reuse.
 3. Please handle and release the cold storage agent gently. If the cold storage agent is broken, please replace it with a new one.
 4. The amount of refrigerating agent in the box/package can be adjusted flexibly to meet the requirements of different customers on temperature and space;
 5. In winter, even if the inner bag and refrigerant are removed, the Portable Refrigerator has a very good thermal insulation effect;
- *Refrigerant: Usually refers to an ice pack, box, or gel ice containing a refrigerant.

BIOBASE GROUP

2# building, No.9 Gangxing Road, High-tech Zone, Jinan City, Shandong Province,
China

Tel: +86-531-81219803/01

Fax: +86-531-81219804

Inquiry: export@biobase.com

Complaints: customer_support@biobase.cc

After-sales service: service_sd@biobase.cc; service_ivd@biobase.cc

Web: www.biobase.cc/www.meihuatrade.com / www.biobase.com